
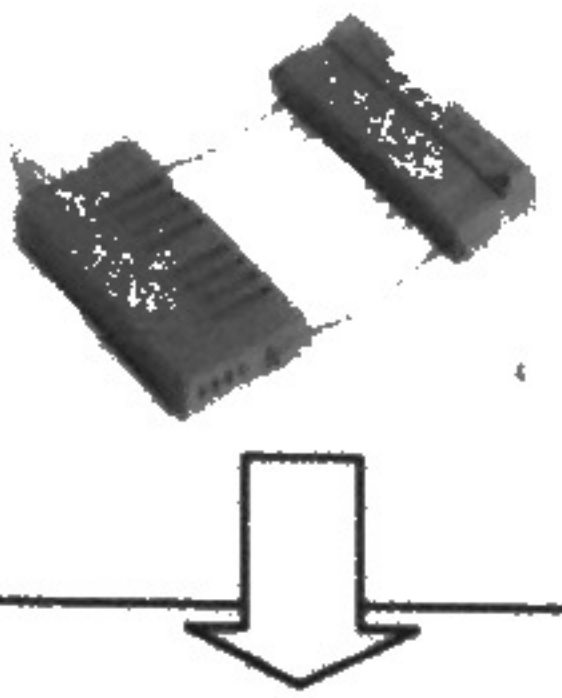
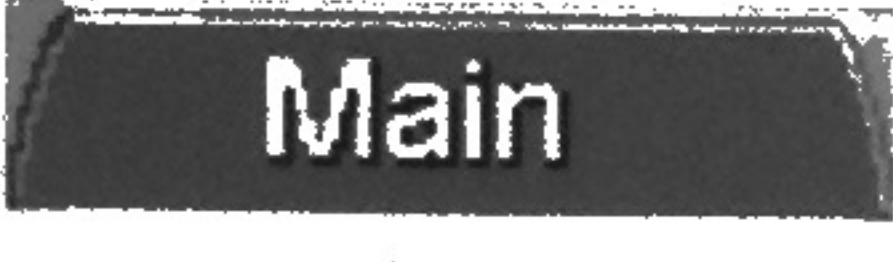
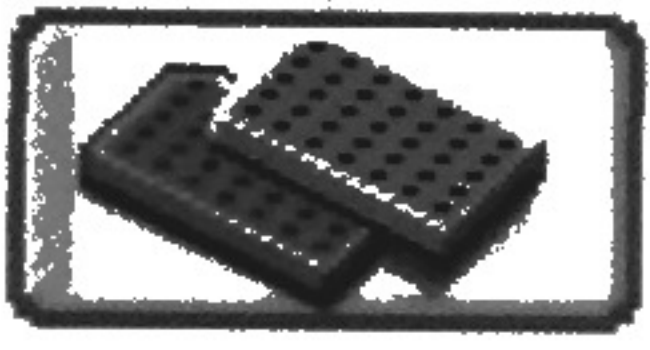
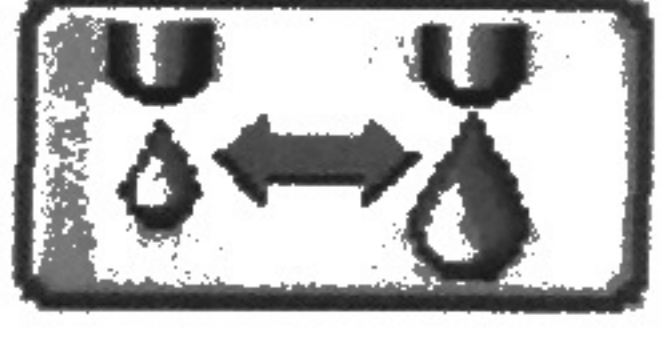



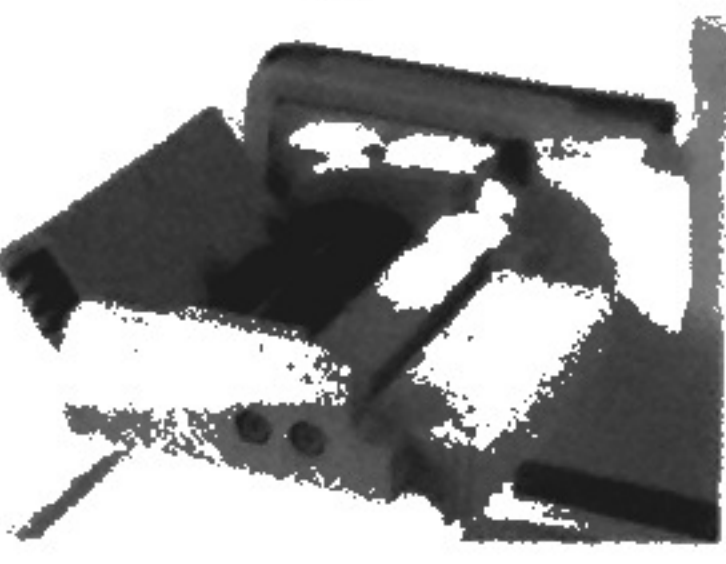




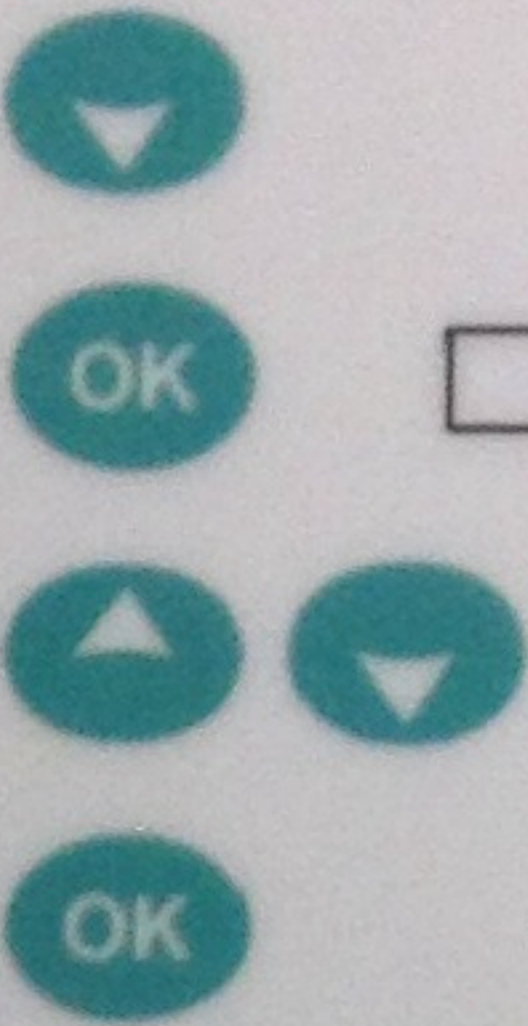


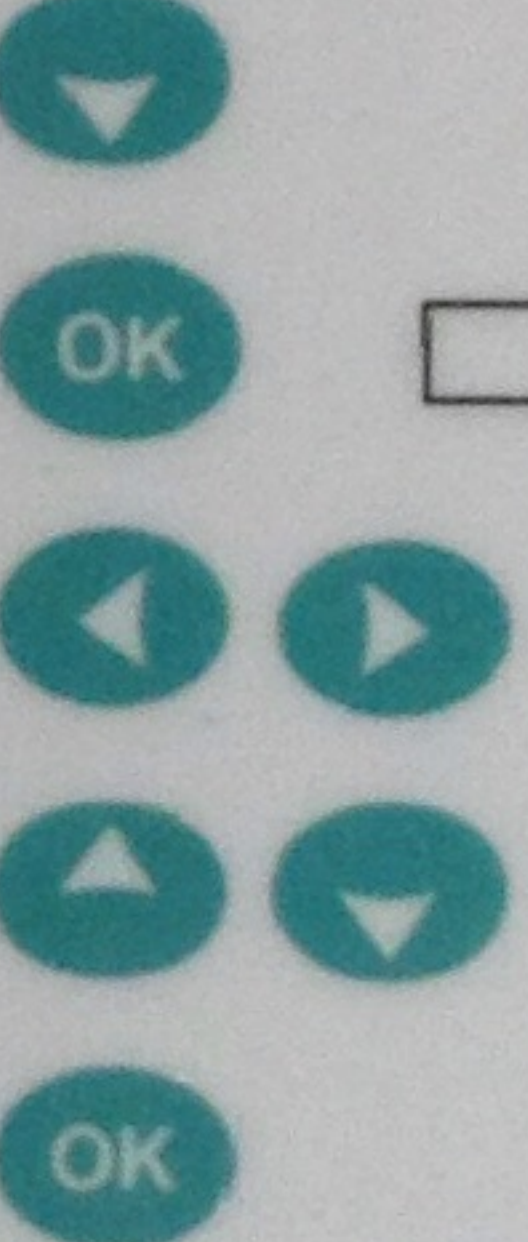
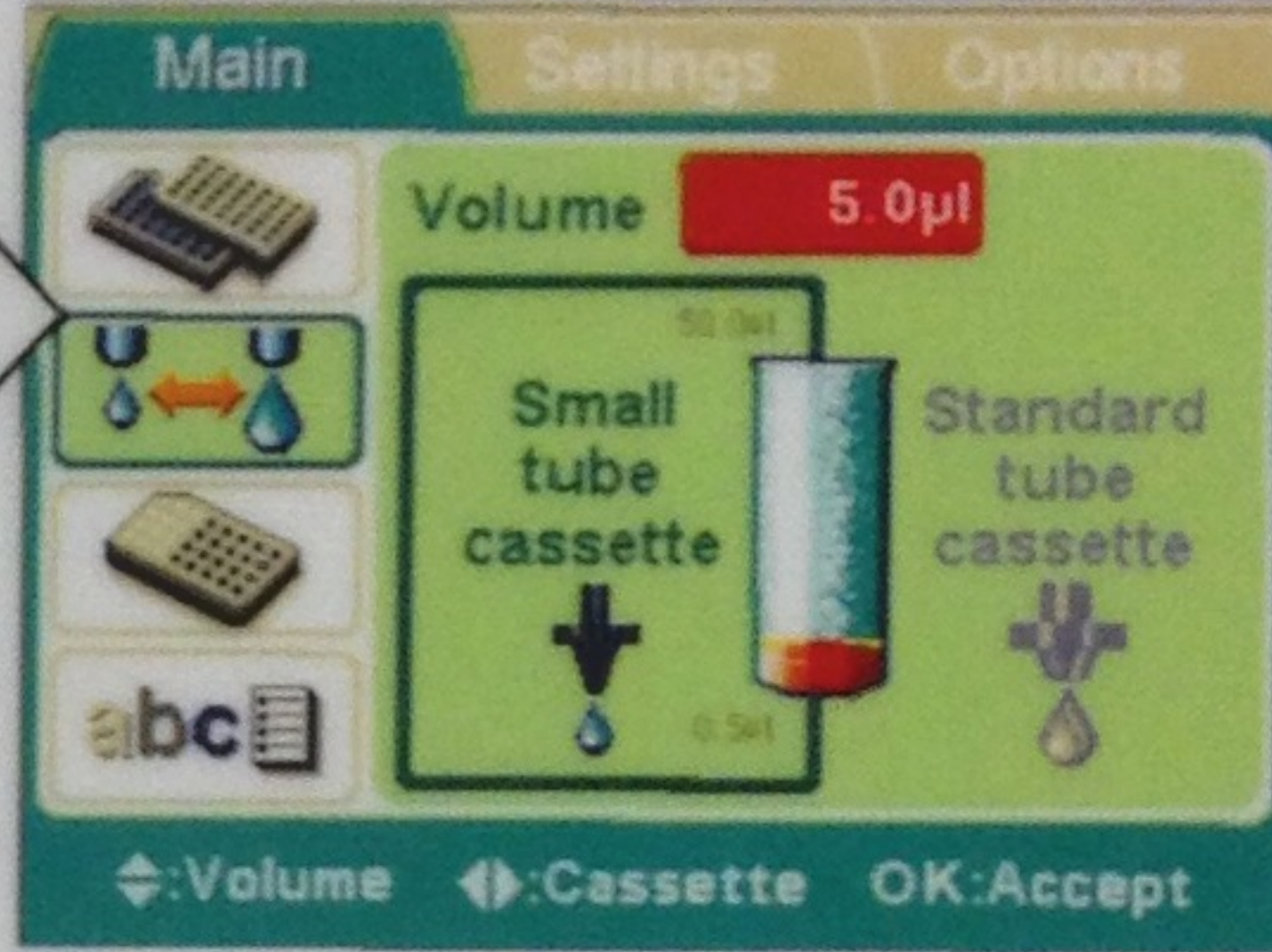
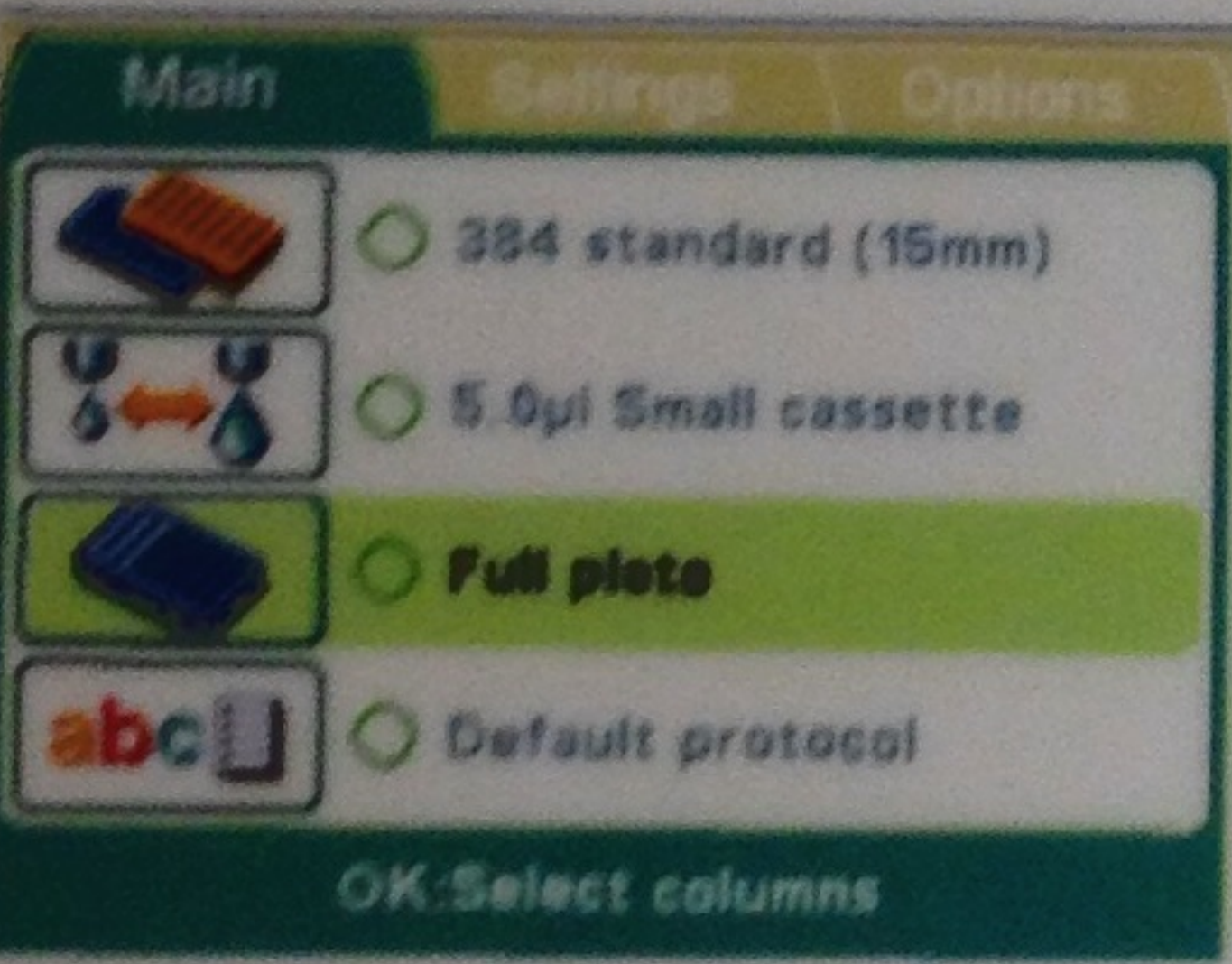


Thermo Scientific Multidrop® Combi

Dispensing Routine dispensing only requires to key in a few buttons.

1.		Power ON the instrument.
2.		Insert the correct dispensing cassette and the priming vessel. Close the rotor cover.
3.		Make sure you are in the Main menu.
4.		Select the plate type and insert the plate.
5.		First select the cassette type and then the dispensing volume.
6.		Keep the PRIME button down until the tubings are completely filled.
7.		Press the START button to start the protocol. The instrument dispenses the plate.
8.		Press the EMPTY button to empty the dispensing tubing.
9.		During prolonged standby, keep the cassette in its rest position. Remove the cassette when the maximum amount of plates are dispensed or you want to change the cassette type in use.
10.		Power OFF the instrument.

Three key selections to start dispensing

After the instrument is turned on, make the following selections to start dispensing when you are in the **Main** menu: plate type, dispensing volume and cassette, and number of columns.

<p>1 - Plate type</p>  <p>OK: Plate type</p>	<p>If the plate type is correct, press OK</p> <p>If you want to change the plate type, press Down Arrow</p> <p>Select the plate type with Left Arrow and Right Arrow</p> <p>To accept, press OK</p>		 <p>⇐: Select OK: Accept</p>
<p>2 - Volume and cassette</p>  <p>OK: Volume and cassette</p>	<p>If the volume is correct, press OK</p> <p>If you want to change the volume, press Down Arrow</p> <p>Select the appropriate cassette with Left Arrow and Right Arrow</p> <p>Set the needed volume with Up Arrow and Down Arrow</p> <p>To accept, press OK</p>		 <p>⇐: Volume ⇐: Cassette OK: Accept</p>
<p>3 - Columns</p>  <p>OK: Select columns</p>	<p>If you intend to fill the whole plate, you can start dispensing after priming the instrument.</p> <p>If you intend to fill only part of the columns, press Left Arrow and Right Arrow</p> <p>To select the columns, navigate with Left Arrow and Right Arrow and to confirm each selection, press OK</p> <p>Last, select Accept with Down Arrow</p> <p>When the Accept box is highlighted, press OK</p> <p>You are now ready to start after filling the tubes through predispense.</p>		 <p>⇐: Select OK: Accept</p> <p>OK: Accept STOP: Escape</p> <p>Selected column(s) = red</p>

Manufactured by:

Thermo Fisher Scientific Oy
 P.O. Box 100, FI-01621 Vantaa, Finland
 Tel. +358-9-329 100, Fax +358-9-3291 0415
 www.thermo.com

Distributed by:

